


# D25<sup>®</sup> NC Packing Instructions

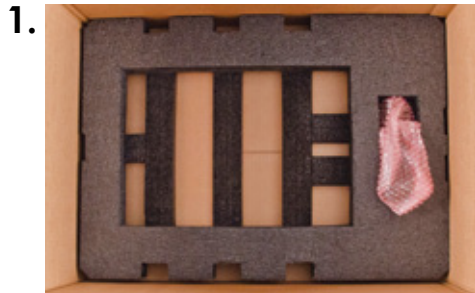

This is the bottom foam section. Place it on the bottom of the box and all the associated cables in the rear pocket of the foam.



Place aluminum pans in center foam cut-out and cover with center foam lid.



Place center foam with lid on top of workstation. Push center foam to back of the sensor until the notch in foam firmly engages with vertical support. Lower sensor tight against center foam lid. Retighten the adjustment handles.




Lift the instrument with center foam.



Place the instrument into bottom foam inside of the shipping box.



Place top foam on sensor. Adjust sensor position as needed to align with cutouts in foam.

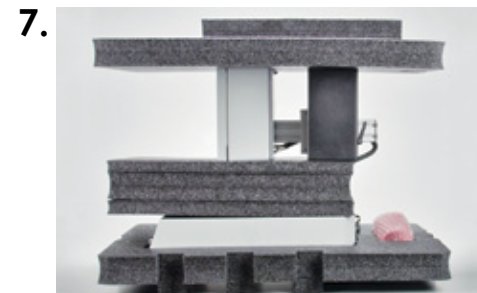

All items are shown as they would sit into shipping box.

Visit [www.hunterlab.com](http://www.hunterlab.com) for information on all our products.