

Event ID	Description	Category
<b>Authentication</b>		
3001	Login(Success): Logged in	Authentication
3002	Login(Failure): Login failed	Authentication
3002	Login failed: Incorrect Account Name and/or Password	Authentication
3002	Login failed: Exceeded the maximum attempts. User Account is locked. Please try to login later.	Authentication
3002	Login failed: The entered user is not a member of Configured Groups.	Authentication
3002	Login failed: Password Expired. Please contact Administrator to change the password.	Authentication
3003	Login attempt: Login tried and cancelled	Authentication
3004	Application logoff: Logged Off	Authentication
3005	Application exit: Exited from Application	Authentication
<b>Sensor Operator</b>		
5001	Add Sensor: Sensor Added	Sensor Operator
5002	Remove Sensor: Sensor Removed	Sensor Operator
5003	Standardization Success: Standardization successful	Sensor Operator
5003	Standardization Cancel/Fail: Standardization Canceled/Failed	Sensor Operator
5004	Sensor Modes: Sensor Mode Added	Sensor Operator
5005	Sensor Modes: Sensor Mode Changed	Sensor Operator
5005	Sensor Modes: Sensor Mode Removed	Sensor Operator
<b>Application</b>		
1001	Modification of Application Preferences: Modified Application Preferences	Application
1002	Modification of Naming Conventions: Modified Naming Conventions	Application
1003	Modification of Default Tolerances: Modified Default Tolerances	Application
1004	Access privileges: Accessed Group Privileges	Application
<b>General</b>		
4003	Read Sample: Sample Read Performed	General
4004	Read Standard: Standard Read Performed	General
4004	Average Selected to New Standard: Standard Read Performed	General
4005	Save Job As: Saved Job File	General
4005	Save Job: Saved Job File	General
4006	Close Job: Closed Job File	General

4009	Sample Record Recall From Database: Sample recalled from Database	General
4009	Standard Record Recall From Database: Standard recalled from Database	General
4009	Series Record Recall From Database: Series recalled from Database	General
4012	Diagnostics test performed: Green Tile Test Performed PASS/FAIL	General
4012	Diagnostics test performed: Repeatability Test Performed PASS/FAIL	General
4012	Diagnostics test performed: Didymium Filter Test Performed PASS/FAIL	General
4014	Universal DB import to Current Job: Imported from Universal DB	General
4014	Universal DB import to New Job: Imported from Universal DB	General
4015	Modified Data Storage: AutoSave Disabled	General
4015	Modified Data Storage: AutoSave Enabled	General
4015	Modified Data Storage: Database Backup Enabled	General
4015	Modified Data Storage: Database Backup Disabled	General
4015	Modified Data Storage: Database Backup Path Changed	General
4015	Modified Data Storage: Database Backup Interval Changed	General
4015	Modified Data Storage: Allow Duplicate Names Enabled	General
4015	Modified Data Storage: Allow Duplicate Names Disabled	General
4015	Modified Data Storage: AutoLogOffTime Interval Changed	General
4015	Modified Data Storage: Database Type Changed to Microsoft SQL Server	General
4015	Modified Data Storage: Database Type Changed to Microsoft Access	General
4015	Modified Data Storage: Database Name Changed	General
4015	Modified Data Storage: Database Server Name Changed	General
4015	Modified Data Storage: Authentication Type Changed to SQL Server	General
4015	Modified Data Storage: SQL Server Username Changed	General
4015	Modified Data Storage: SQL Server Password Changed	General

