



Reflectance Sample Shelf with Light Cover (B02-1005-172)

Instrument: ColorQuest XE, UltraScan PRO, UltraScan VIS

Purpose: The optional reflectance sample shelf with light cover can be placed at the reflectance port for measurement of powders, granules, pastes, or liquids held in a Transmittance cell (10-mm, 20-mm, 33-mm, or 50-mm cell). An opaque cover is also available to prevent ambient room light from affecting the measurement.

Installation:


1. Lower the sample clamp as far as it will go.
2. Insert the two finger clamps of the sample shelf into the guidance holes located below the reflectance port on the front of the instrument.



3. Tighten the sample shelf into place by turning both knobs clockwise.

Sample Measurement:

1. Place your sample on the shelf and cover it with the light cover (if included).

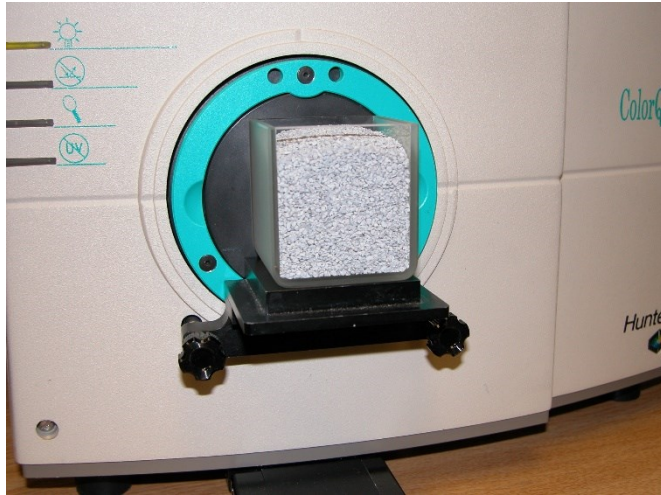

2. Measure the sample.