



4-mm Port Plate (A02-1011-184)

Instrument: ColorQuest XE, UltraScan PRO, UltraScan VIS

Purpose: This option provides a modified port plate with a 4-mm round opening. This port plate should only be used when the small area view option is available and in use.



Installation: To install the port plate, remove the standard port plate and snap the modified port plate into place at the reflectance port.

Sample Measurements: Standardize the instrument in a reflectance mode using the 0.375-inch (9.525 mm) area view.

Note: *This port plate has an opening that is smaller than the usual 1-inch port plate. Performance specifications (i.e., factory-supplied color values, stability, accuracy) are based on the standard 1-inch round port plate. These specifications will not be met with the modified port plate. The change of geometry that occurs will cause a slight shift in the color values of the calibration standards.*