



Glass Port Insert

Instruments: ColorFlex EZ (02-6624-05), ColorQuest XE (A01-2022-124), LabScan (A02-1010-319), UltraScan PRO & UltraScan VIS (A01-2022-124)

Purpose: This option provides a snap-in port plate with an opening one inch in diameter that is covered with glass. The glass insert protects the inside of the instrument from sample and airborne particles.

Installation: Remove the standard port insert by pulling it out of the port insert retainer. Replace it with the glass port insert. Make sure that the flat portion of the insert is facing outward with the beveled portion toward the interior of the sensor.

Sample Measurements: Standardize the instrument with the glass in place. Check the measurements of the green and white tiles before performing any other diagnostics without the glass cover. Difference measurements using a glass cover will be more accurate than absolute measurements.

