



## Sample Clamp and Port Forward Stand

(D2-1010-459)

**Instruments:** ColorFlex EZ

**Purpose:** The port-forward stand holds the instrument securely in the port-forward position and the sample clamp holds samples flat at the sample port.

**Installation:** To attach the stand to the ColorFlex EZ, lay the instrument on its side and install the two screws through the stand into the bottom of the instrument using the Allen wrench provided.

**Sample Measurement:** To use the sample clamp, pull the spring-loaded clamp away from the sample port, insert the sample, and gently release the clamp.

