



Skein/Swatch Holder Option

(02-7396-00)

Instruments: ColorFlex EZ, LabScan, MiniScan EZ, ColorQuest XE, UltraScan Pro, UltraScan VIS

Purpose: This is a device for measuring yarn skeins. Wind the yarn around the skein holder in multiple taut layers until it is effectively opaque and is as flat as possible. Secure it in place with the detachable arms on the sides of the skein holder.

Sample Measurement: Place the skein holder at the sample port and back it with the sample clamp or a white backing tile to provide a consistent background and pressure. Make several measurements of the skein, rotating the holder 90° between measurements and averaging the readings for the final result.

