



45/0 LAV Nose Cone with Lower Glass Assembly (A02-1002-129) or Nose Cone with Lower Polycarbonate Assembly (B02-1009-500)

Instruments: MiniScan

Purpose: This option includes a special nose cone with a removable glass or plastic port cover assembly. The nose cone (black) portion is intended for permanent use on the instrument and should not be removed. Replacement cover assemblies (the lower portion of the device) are available from HunterLab. The assembly is sealed, but is not to be considered waterproof.



Glass Nose Cone



Replacement Glass Cover



Nose Cone Installed

Installation: The nose cone/port cover is generally in place when the MiniScan EZ is shipped from the factory. However, if installation is required, place the assembly over the instrument port and secure it using three Phillips-head screws. If the glass or plastic cover is to be replaced, remove the three machine screws with lock washers, replace the window and the O-ring, and secure it using the machine screws.