



UV Port Insert

(02-6624-05)

Instruments: ColorFlex EZ

Purpose: This port insert contains a 420-nm filter to prevent light below 420 nm from hitting the sample. This is important when samples have been treated with UV enhancing chemicals.

Installation: Remove the standard port insert by pulling it out of the port insert retainer. Replace it with the UV port insert. Make sure that the flat portion of the insert is facing outward with the beveled portion toward the interior of the sensor.

Sample Measurement: Standardize the instrument with the UV filter in place. Check the measurements of the green and white tiles before performing any other diagnostics without the UV filter.