



Tomato Red Tile

(L02-1014-594)

Instruments: ColorFlex EZ

Purpose: This option provides an uncalibrated tomato red porcelain enamel on steel instrument tile. Tomato score values may be assigned to the tile based on known USDA tomato color standards to hitch the instrument to those soft standards. Alternately, the TILE-CAL-TRACEABLE option can be purchased from HunterLab to have tomato score and color values assigned to the tile at the factory.

Sample Measurement: When a tomato score is to be used, read the tomato tile as a standard and hitch the instrument to the L, a, b values given on the standard label. Save the standard and recall it to re-hitch the instrument whenever tomato scores are to be measured.