



420-nm UV Filter Assembly (A02-1009-954)

Instrument: MiniScan

Purpose: This option provides a 420-nm UV filter in a replaceable assembly for 45/0 LAV MiniScan EZs. The UV filter may be replaced or removed when necessary; however, the instrument is not to be used without the special nose cone in place.

Note: *The special nose cone required for use of this part is not included, but may be purchased separately.*

Installation: To install the UV filter assembly, secure it to the instrument nose cone using three 4-40 pan-head screws with split-lock washers.

Replacement UV filter assemblies are available from HunterLab and are marked “UV” to differentiate them from cover glass assemblies. To replace the UV filter assembly, remove the screws and lift the assembly off the nose cone. Place the new UV filter assembly on the nose cone and replace the screws.