

# Addendum to ColorFlex EZ User's Manual

---

## Description

CMR-2570 provides a fiber package adapter for use when measuring bobbins using the ColorFlex EZ instrument. The adapter should be ordered based on the diameter of the bobbin to be measured, referencing the following part numbers:

Part Number	Bobbin Diameter (inches)	Bobbin Diameter (mm)
CMR-2570-P1	6	152
CMR-2570-P2	7.5	191
CMR-2570-P3	9-10	229-254
CMR-2570-P4	5	127
CMR-2570-P5	12-14	305-356
CMR-2570-P6	3.5	89



## Installation

1. To install a fiber package adapter, first remove the standard ColorFlex EZ port insert. Note the alignment hold on the display side of the measurement port.
2. With the alignment pin on the port adapter directly above the alignment hole near the instrument port, snap the port adapter desired into place over the port.

## Operation

Standardize the instrument with the standard port plate in place and then install the fiber package adapter as described above. To measure a bobbin, simply place the bobbin on the port adapter. You may need to hold it in place. Averaging several measurements of different areas of the bobbin is recommended.

