


## Addendum to UltraScan PRO User's Manual

---

### Description

CMR 3030 provides a small area view (7-mm/0.25-inch port size) port plate with cover glass.

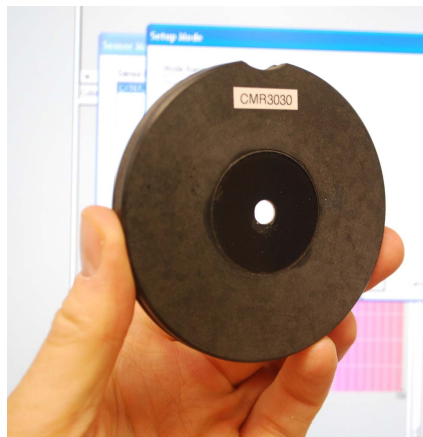

### Installation

Remove the existing port plate and replace it with the CMR 3030 port plate.

### Operation

Standardize the instrument for the 0.190-inch (4.826-mm) area view with the CMR 3030 port plate in place. Then proceed with sample measurements using the CMR 3030 port plate.

*Note: Absolute values read with the glass port plate in place may not be identical to those read with the regular port plate.*