

Addendum to ColorQuest XE SAV, UltraScan VIS and UltraScan PRO User's Manual

Description

CMR-3110 provides an additional holder for the D02-1012-931 Semi-Micro Powder Holder used for the color measurement of small quantities of packed powder. The additional CMR-3110 parts allow multiple samples to be prepared in advance.

Installation/Operation

1. Place the powder to be read into the powder holder.
2. Pack the powder down using the plunger. Do not overtighten the powder holder. The powder should be at least 1/4-inch deep in the holder.

