

Addendum to User's Manual

Description

This CMR modification incorporates a spill tray with a standard C02-1005-481 transmission cell holder and an L-shaped bracket that centers a disposable, plastic 10-mm transmission cell at the TTRAN port of any of HunterLab's sphere instruments— ColorQuest XT, ColorQuest XE, UltraScan VIS and UltraScan PRO.



Installation

Align the holder against the TTRAN port on the instrument and tighten the screws to secure the device to the transmission floor.

Operation

This device is designed to fit tightly at the TTRAN port of all HunterLab sphere instruments minimizing potential damage should a spill occur.

