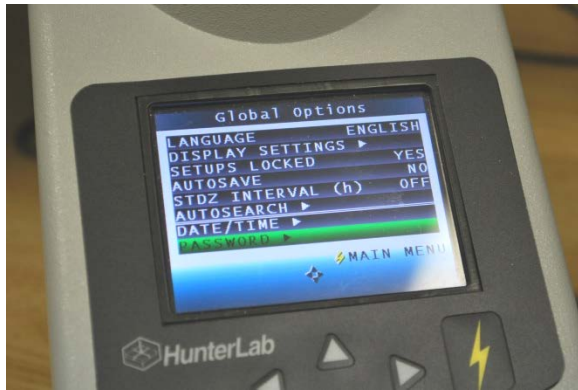


Addendum to ColorFlex EZ User's Manual

Description

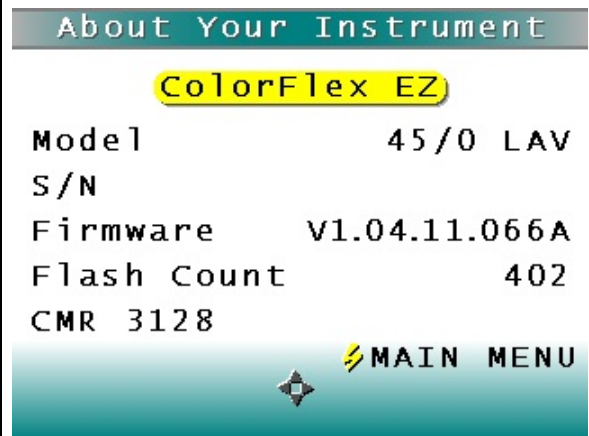


This CMR modification to the ColorFlex EZ firmware adds password protection before product setups may be modified, reset or recalled. This option is typically installed at factory.

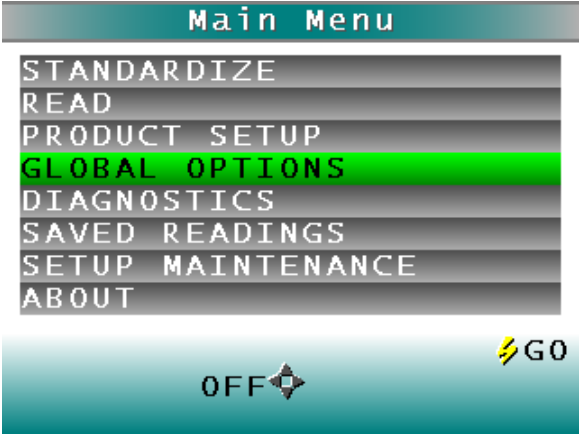

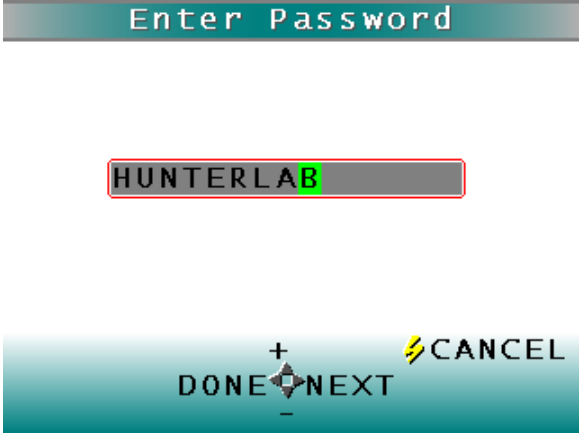
Installation

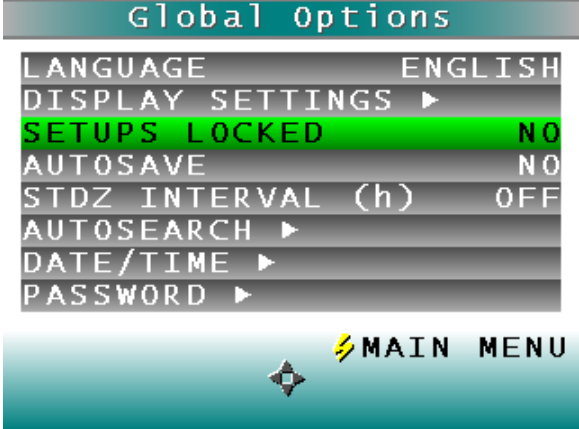
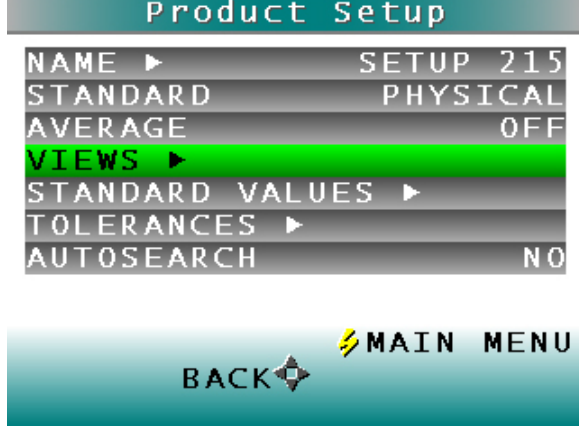
The special CMR-3128 firmware for the CFEZ will typically be installed at the factory.

Go to MAIN MENU/HELP/ABOUT and you should see CMR 3128 listed.



Operation

<p>When CMR3128 is factory installed, the instrument will typically ship with the Main Menu/Global Options/Setups Locked = “YES” and the default Password set to “HUNTERLAB”. Before any setups can be modified, the User should change the password to one of their choosing.</p>	
<p>To reset the password, go to MAIN MENU/GLOBAL OPTIONS/PASSWORD.</p>	
<p>Enter the password “HUNTERLAB” to allow access. Then enter a new password of your choice.</p> <p>Note: Spaces at the beginning or end of an entered password are ignored.</p>	

<p>To unlock the setups, go to MAIN MENU/GLOBAL OPTIONS/SETUPS LOCKED and select change “YES” to “NO”. You will be prompted for your new password to make this change.</p>	 <p>The screenshot shows the 'Global Options' menu. The 'SETUPS LOCKED' option is highlighted in green and set to 'NO'. Other options include LANGUAGE (ENGLISH), DISPLAY SETTINGS, AUTOSAVE (NO), STDZ INTERVAL (h) (OFF), AUTOSEARCH, DATE/TIME, and PASSWORD. A 'MAIN MENU' button is at the bottom.</p>
<p>From the Main Menu, select a Product Setup to modify.</p>	 <p>The screenshot shows the 'Main Menu' with 'PRODUCT SETUP' highlighted in green. Other options include STANDARDIZE, READ, GLOBAL OPTIONS, DIAGNOSTICS, SAVED READINGS, SETUP MAINTENANCE, and ABOUT. A 'GO' button is at the bottom right.</p>
<p>Within a Product Setup, enable a View and make any required changes to the setup.</p>	 <p>The screenshot shows the 'Product Setup' menu. The 'VIEWS' option is highlighted in green. Other options include NAME (SETUP 215), STANDARD (PHYSICAL), AVERAGE (OFF), STANDARD VALUES, TOLERANCES, and AUTOSEARCH (NO). A 'MAIN MENU' button is at the bottom right.</p>

To lock the setups, go to MAIN MENU/GLOBAL OPTIONS/SETUPS LOCKED and select change “NO” to “YES”. You will not be prompted for a password to lock the setups; only to unlock.



Whenever the instrument is turned on or comes out of stand-by mode (instrument has power but screen is dark), the setups revert to locked status (GLOBAL OPTIONS/SETUPS LOCKED = “YES”).

Unless setups are unlocked by password, new setups cannot be uploaded into the sensor from a flash disk or EasyMatch QC software or the setups reset under GLOBAL OPTIONS/SETUP MAINTENANCE.

If a User forgets their installed password, a backdoor option is available by contacting HunterLab at support@hunterlab.com.