

Addendum to ColorFlex EZ User's Manual

Description

CMR-3200 provides a 63mm diameter glass round that fits in the ColorFlex EZ Open Port Insert for 2.5 in (64 mm) Sample Cup (HunterLab part number 04-6622-00). This glass round is used during standardization to negate the absorbance effects of the sample cup window.

Installation

1. Install the ColorFlex EZ Open Port Insert for 2.5 in (64 mm) Sample Cup (HunterLab part number 04-6622-00).
2. Place CMR-3200 glass round on the port insert.

Operation

1. Install the ColorFlex EZ Open Port Insert for 2.5 in (64 mm) Sample Cup (HunterLab part number 04-6622-00).
2. Place CMR-3200 glass round on the port insert.
3. Standardize the instrument as described in the user's manual.
4. Remove CMR-3200 glass round and proceed with sample measurements using the glass sample cup (HunterLab part number 04-7209-00).

