



ColorFlex EZ Sensor Port-Down Stand CMR-3220

Instrument: ColorFlex EZ

Purpose: The sensor stand permits port-down mounting of the sensor.

Installation: To install the instrument onto the stand, turn it with the port down and then insert the knob screws through the back of the stand and into the receiving holes on the back of the ColorFlex EZ.



Figure 1. Stand with Knobs



Figure 2. ColorFlex EZ Ready to Read

The unit can be raised and lowered, and for a series of samples, the distance between the surface of the sample and the instrument port need to be zero, that means that the sample needs to be in contact with the instrument port.

Using EasyMatch QC software, the instrument can be standardized from this position and samples can be placed into a container below the instrument port for reading.

Note: *The sample in the container should be in contact with the port.*



Figure 3. Port-Down Sample Measurement